

# 012656 SEC-2F/3F Au grid electrode for flow cell

---

Thank you for purchasing an Au grid electrode. Please read the following instructions carefully before using the Au grid electrode.

We do not guarantee any use against these precautions.

1. Do not sonicate the Au grid electrode.
2. Do not polish the Au grid electrode.
3. Do not use the Au grid electrode in a strong acid or a strong alkaline solution, and do not clean it with those solution.
4. Do not rub the surface of the Au grid electrode with a hard material, or it may cause the disconnection of the electrode part.
5. Do not use the Au grid electrode under a higher or lower temperature.
6. Any modification on the surface of the Au grid electrode with chemicals or physical treatments is the responsibility of the customer. The characteristics of the modified electrode cannot be guaranteed.
7. Do not apply a higher or lower potential to the Au grid electrode.
8. The Au grid electrode is made on the glass, so it may be damaged by unreasonable force such as twisting.
9. Do not touch the surface of the Au grid electrode with fingers, or a grid pattern may come off or break.
10. Do not clean the Au grid electrode with an ozone cleaner.

## CAUTION:

Except with the serious defect of Au grid electrode due to our production process, the electrode can not be replaced with a new one. In addition to that, minor not serious matter such as pinholes in the lead and glass burrs is not the subjects of warranty.