

012657 SEC-2F/3F Carbon grid electrode for flow cell

Thank you for purchasing a Carbon grid electrode. Please read the following instructions carefully before using the Carbon grid electrode. We do not guarantee any use against these precautions.

1. Do not sonicate the Carbon grid electrode.
2. Do not polish the Carbon grid electrode.
3. Do not use the Carbon grid electrode in a strong acid or a strong alkaline solution, and do not clean it with those solution.
4. Do not rub the surface of the Carbon grid electrode with a hard material, or it may cause the disconnection of the electrode part.
5. Do not use the Carbon grid electrode under a higher or lower temperature.
6. Any modification on the surface of the Carbon grid electrode with chemicals or physical treatments is the responsibility of the customer. The characteristics of the modified electrode cannot be guaranteed.
7. Do not apply a higher or lower potential to the Carbon grid electrode.
8. The Carbon grid electrode is made on the glass, so it may be damaged by unreasonable force such as twisting.
9. Do not touch the surface of the Carbon grid electrode with fingers, or a grid pattern may come off or break.
10. Do not clean the Carbon grid electrode with an ozone cleaner.

CAUTION:

Except with the serious defect of Carbon grid electrode due to our production process, the electrode can not be replaced with a new one. In addition to that, minor not serious matter such as pinholes in the lead and glass burrs is not the subjects of warranty.