

013338 DRE-GCD GC disk

Thank you for your purchase from BAS Inc.

Please read the following precaution before using the products.

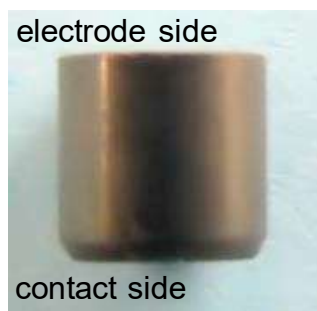


Fig.1 Electrode direction



Fig.2 Label of serial code

This product can be used with 4 mm diameter disc electrode of DRE disc replaceable electrode.

Pay attention the direction of disk electrode at assembly(Fig.1).

The chamfered side is the back "contact side".

You can check the assembly, disassembly and polishing method of DRE disk replaceable electrode on our website.

<https://www.als-japan.com/1571.html#defaultTab15>

013608 DRE-DCP Disk electrode polishing and exchanging tool kit must be used for assembly, disassembly or polishing of DRE electrode.

Damage caused by using other tools is not covered by our warranty.

CAUTION

1. Be sure to keep the electrode label(Fig.2). The serial code is required for any case of inquiry or confirming the checking data of electrode.
2. Do not soak DRE disk replaceable electrode in strong acid or base solution for long time. After use, disassemble and wash electrode with pure water immediately.
3. Use the electrode at room temperature and atmospheric pressure.
4. The electrode should be protected from strong shock.
5. Do not soak the electrode in organic solvent for a long time.
6. We cannot replace the product after using.
7. Do not use damaged electrode.
8. The disk modification by chemical or physical process will change the electrode surface state. In this case, the electrode surface regeneration is not warranted.

You can browse the checking data of electrode in below URL .

<https://www.als-japan.com/dl/>

Please check color pdf data and the latest information with our website.

BAS Inc.

<https://www.als-japan.com> email: sales@als-japan.com