



SEC2000 Ver1.3 Quick-Start Guide

Shipment Components

The following information and documentation ship with the SEC2000 Spectrometer:

1. Software and Manual Resources USB memory

The following software and manual are included in the USB memory.

- ① Visual Spectra 2.1 Sr (Software installer, Driver and Manual update installer for WinXP/Win7)
- ② SEC2000 Manual (Operation manual)
- ③ SEC2000 Ver1.3 Quick-Start Guide

2. Wavelength Calibration Data Sheet

3. USB cable

4. Trigger cable

5. Power supply for light source

6. SEC2000-DH-RNG (only with system)

Installation Sequence is Important for Success

Setup is easiest if you use the sequence below.

1. Install the software Visual Spectra 2.1 and Device Driver on PC.

Insert the USB memory to your computer and install the software by using the setup program. Then, install the Driver.

2. Connect the SEC2000 spectrometer to the PC by USB cable.

Using the enclosed USB cable, connect the spectrometer to the PC. The PC automatically recognizes the USB interface when you use Win XP. Please follow the steps displayed.

When the tasks have been completed, the SEC2000 spectrometer is ready to use.

You can run experiments and gather data with your Visual Spectra 2.1 software.